- Zum Nachfüllen nur Original Kigas-Kartuschen von Kisag verwenden.
- N'utilisez qu'une cartouche de gaz originale Kigas pour le remplissage.
- Per il riempimento, utilizzate esclusivamente delle bombole originali Kigas.
- For refill use only original Kigas canisters of Kisag.

Art. 8400

Art. 8401

Kisag AG
 CH-4512 Bellach
 Tel. +41 32 617 32 60
 Fax +41 32 617 32 70
 E-Mail: kisag@kisag.ch
www.kisag.ch

3.2013/2800.902

DE

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EN

Kisag Bunsen Burner
 Operating Instructions

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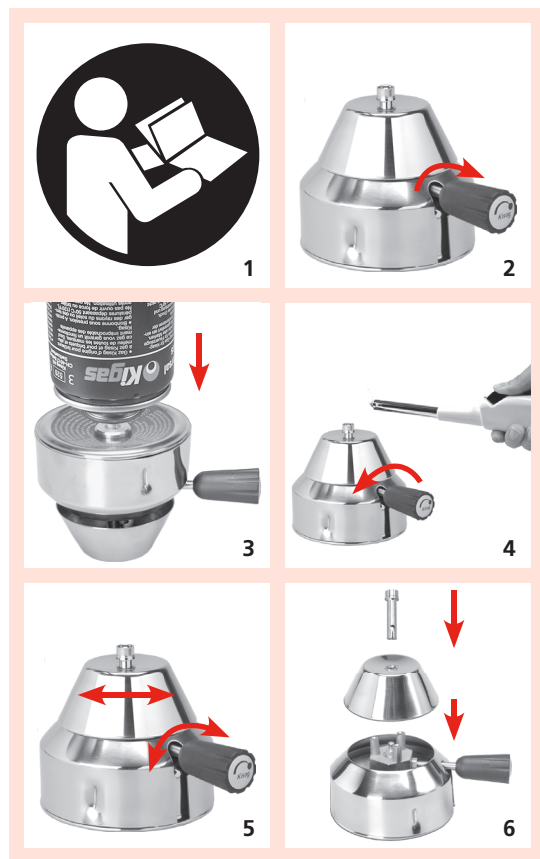
EM | USE ACCORDING TO THE PURPOSE

The Bunsen burner of Kisag, type 9500 is a device of one flame and is designed for use in laboratories. The conditions for safe operation are to comply with all safety rules and operating according to this manual.

Manufacturer	Kisag AG, 4512 Bellach, Switzerland
Kind of gas	Butane/Propane mixture
Kigas refill cartridges 400 ml	Lasts for more than 6 fillings
Category of appliance	Direct pressure of butane/propane gas mixture
Nominal thermal load	70 g/h; approx. 1 kW
Burning duration	Approx. 2 h/filling (working flame)
Temperature	Until 1300° C
Jet	No. 21

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EM | GUARANTEE

Each damage on the appliance caused by faulty materials or faulty processing, which appears within two years after date of the purchase document, will be replaced free of charge. For damages caused by careless use, use of defective parts, repairs by third parties, normal wear and tear or for any consequential damages there is no claim for guarantee or responsibility existing. Immediately after statement of a damage the appliance must be sent together with a copy of the purchase document and a description of the damage stated, transport fees paid, to Kisag AG, CH-4512 Bellach, or to an after-sales service licensed by Kisag.